



# Lipid Staining Kit

Cat# ST-R100

## INSTRUCTION MANUAL (ZBM-12)

### STORAGE CONDITIONS

- **Reagents & Buffers:** Room temperature 25°C

**For *in vitro* Use Only**

***For Research Use Only. Not For Use In Diagnostic Procedures***

### LIMITED PRODUCT WARRANTY

This warranty limits our liability to replacement of this product. No other warranties of any kind, express or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Zen-Bio, Inc. Zen-Bio, Inc. shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

### ORDERING INFORMATION AND TECHNICAL SERVICES

- **Zen-Bio, Inc.**
- **3200 Chapel Hill-Nelson Blvd., Suite 104**
- **PO Box 13888**
- **Research Triangle Park, NC 27709**
- **Telephone** (919) 547-0692
- **Facsimile (FAX)** (919) 547-0693
- **Toll Free** 1-866-ADIPOSE (866)-234-7673
- **Electronic mail (e-mail)** [information@zen-bio.com](mailto:information@zen-bio.com)
- **World Wide Web** <http://www.zen-bio.com>

## Items included in the kit:

The protocol is designed for assay of cells in a 96-well format. For other formats, please adjust the volumes added to each well according to the surface area of the well/flask you are using.

ITEM	STORAGE NOTES	UNIT	QTY
Fixative Solution	Store at room temperature, preferably in a cabinet approved for storage of flammable chemicals.	40ML BOTTLE	1
Oil Red O Stock Solution	Store at room temperature in a dark bottle, preferably in a cabinet approved for storage of flammable chemicals.	6ML BOTTLE	1
Syringe	10 cc syringe	EACH	1
Filter	1µm PTFE filter	EACH	1
Tray	For multi-channel pipetters, clear polyvinyl	EACH	3

### Other equipment/reagents required but not provided with the kit:

- Mature adipocytes or other lipid-containing cells to be stained
- Deionized or distilled water (dH<sub>2</sub>O)
- Multi-channel Pipet , single channel pipet and pipet tips

## Assay Procedure

### 96-well plate format Protocol

1. Remove row by row all the liquid and add 100µl/ well Fixative Solution. Allow the plate to set at room temperature for about 5 minutes. Repeat liquid exchange with another 100µl/ well Fixative Solution.
2. Cells must be fixed for at least 1 hour at 4°C. Alternately, cells in Fixative Solution may be sealed using sealing film/foil and kept at 4°C for at least 7 days before proceeding.
3. **Prepare Working Oil Red Solution (WOROS): (Prepared fresh on day of assay) Prepare a working solution 40% water and 60% Oil Red O stock. Mix by inversion.**
  - a. This working solution must be kept at room temperature at least 20 minutes before filtering.
  - b. Filter out particulate material using a 1µm polypropylene PTFE syringe filter (supplied with the kit) or #1 Whatman filter paper. NOTE: Filter paper is effective; it just takes longer to filter. Polycarbonate filters will dissolve in isopropanol; DO NOT USE THEM.
4. Remove all the Fixative Solution. Wash at least twice using a multichannel pipetter by adding 200 µl/well dH<sub>2</sub>O. Remove all the water using a multichannel pipetter. Make sure all the water is removed BEFORE proceeding.
5. Add 50 µl/ well Working Oil Red O solution (WOROS) at room temperature for 10-15 minutes. Be careful not to touch the sides of the wells; the pipet tips should go at a 90° angle into the wells.
6. Remove all the Oil Red O solution from each well. Wash 2-3 times with 60-80µl/ well dH<sub>2</sub>O. Pictures may be taken at this time. Leave water inside wells during photography.